





Atom Physics

Instadose App Setup



1. Open your preferred app store on your mobile device and search for the 'Instadose' app. Select the download button.
2. Your device's Bluetooth must be turned on to transfer data through the app.
3. Open the app and enter your username and password.

Note: If you forget your password, you can reset it or contact Atom Physics.

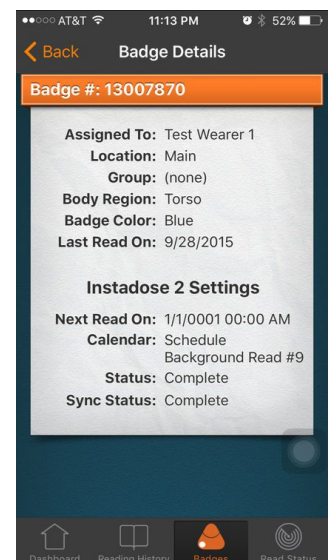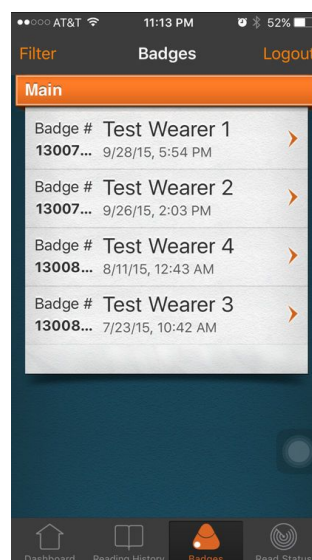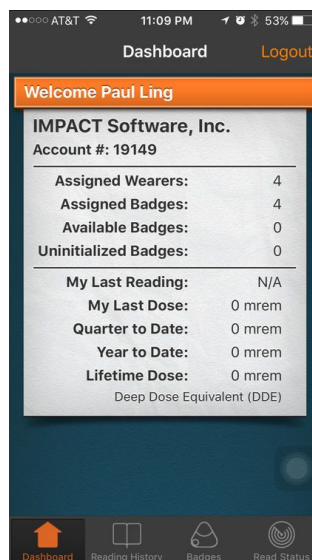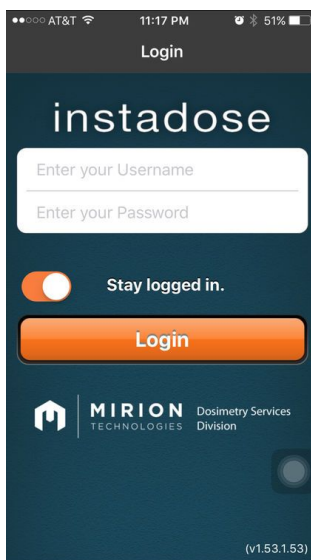

The Dashboard, Badges, and Badge Details screens can be used to review your active badges.

For help with the Instadose app, contact Atom Physics at 303.748.5499 or info@atomphysics.com

Atom Physics

Instadose Badge Reading



Your badge has already been initialized and is constantly collecting data. However, logging readings in the app is the only way to see the dose exposure readings collected by the badge. It is recommended to do this monthly.

1. Open the instadose app (with Bluetooth on) and select the 'Read Status' icon on the bottom of the screen
2. Hold the badge near your device. Press and hold the button on the back of the badge for 5 seconds until the indicator light stays green for one second.
[Note: If a red light appears, wait one minute then try again. If the problem persists, contact Atom Physics at 303.748.5499 or info@atomphysics.com for troubleshooting.](#)
3. Wait up to 60 seconds for all of the dose readings to transfer to your device. The indicator light will stay green for one second when it is complete.
[Note: The read status screen on your device will not confirm the data transfer.](#)
4. To verify a successful communication, click on the 'Read History' icon on the bottom of the Dashboard. You may need to refresh the history (e.g. pulling down on the iPhone). Make sure recent dose readings appear.
[Note: An asterisk \(*\) denotes a reading lower than 1 mrem](#)

